CWE 7745 /CWB 7726
Bedienungsanleitung
Operating Instructions
Mode d'emploi
Modo de utilização
Manual de uso

CE

SOEHNLE

PROFESSIONAL

Ihr Soehnle-Servicepartner

Your Soehnle-Partner

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Technical Information

- Operated by means of a foil keypad.
- Mains operation with power supply unit (Mod. 7745) or incorporated power supply (Mod. 7726)
230 V/50-60 Hz/ca. 12 VA
(battery-/accu-operation see instructions).
- Operating temperature – 10° +40° C.
- Storage temperature – 40° +70° C.
- Manual resetting range – 1 % +3 %.
- Initial loading possible (Please contact your dealer).
- All data is stored securely against power failure.
- Serial RS 232 (V24) interface.

Specification of the nameplate

	<p>I↔I1 Weighing range (weighing ranges in the case of multi-range scales)</p> <p>Max Range permitted for calibrated</p> <p>Min weighing</p> <p>e= verification scale internal</p> <p>SNR Serial number (scales type, last digits of year of manufacture, number)</p> <p>CEXX EU-conformity symbol indicating year of attachment, corresponding to the beginning of verification validity (e.g. CE98).</p> <p>Ⓜ accuracyclass, EC Pattern approval (e.g. D98-09-002).</p> <p>XXXX Code of the office quoted (this office stamped the scale initially).</p> <p>M EC metrology mark</p>
--	--

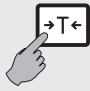
Getting started

	<p>Set up the weighing platform at the place required. Use the water level and the adjustable feet to ensure that the scale is positioned horizontally.</p>
<p>kg lb</p>	<p>Kg/lb-conversion</p> <p>Press key first and hold, then press key additionally.</p>

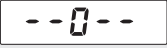
Display symbols

	Zero message
	Idle indicator
	Active range in the case of multi-range scales
NET	Net weighing
	Battery operation: Replace battery respectively charge accu.

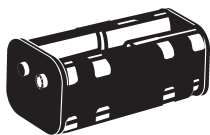
Using the scale

  	<p>Weighing</p> <p>Switch on (only when the scale is not loaded).</p> <p>After the test routine, the main display is set to zero.</p> <p>The scale is ready for use.</p> <p>Place the items on the scale.</p>
 	<p>Manual zero resetting</p> <p>When display does not show zero (scale unloaded). Manual zero resetting range see "Technical information".</p>
 	<p>Weighing with tare</p> <p>Determine the tare of the empty container. Scales are ready to perform net weighing.</p>
	<p>Deleting the tare weight</p> <p>The gross weight appears in the display or "0" when the scale is unloaded.</p>
	<p>Printer</p> <p>Press key to print. Your service partner can modify the printout.</p>
 	<p>Switch off</p> <p>Possible immediately when display shows "0". During weight indication keep pressing for 3 seconds.</p>

Malfunctions

	Indicator	Remedy
	An acoustic signal confirms that entry was correct. Incorrect entries are indicated by 5 acoustic signals in swift succession.	
	The scale sets the zero value automatically when it is switched on. If the scale is outside the tolerance range --0-- is.	Remove the items from the weighing platform. Remove any dirt. Check whether weighing platform is mounted duly.
	Underload: Only the bottom horizontal lines appear on the display.	Switch the scale off and on again. The zero point will be reset automatically.
	Overload: Only the top horizontal lines appear in the display.	Remove some of the items to be weighed from the scale.
   	The zero setting range has been exceeded. Taring is not possible in case of over- and not loaded balance. Printing is not possible in case of over- or underloaded balance. Conversion to unit lb not admissible.	
		In case of all other error indications: Please contact your service partner.

Accu-/battery operation

At the bottom of the scale you can find the battery holder. 8 batteries "baby cells 1,5 C" or NiCd-accumulators of the same size are required. The length of operation depends on the quality of the batteries or accumulators, e.g. at an accu-capacity of 1,4Ah appr. 8-10 operation hours. During accu-resp. battery operation the background light of the display is switched off automatically.

If rechargeable batteries are not used for a longer time, they must be recharged every 4-5 months.

If accumulators are used they cannot be recharged in the scale.

Battery chargers are available in electronic stores.

	Message during battery operation	Remedy
	Residual length of operation appr. 10 min.	
	Display shows Battery-empty-message.	Change batteries or recharge accumulators.
	Scale cannot be switched on during accu or battery operation.	Change batteries or recharge accumulators.

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SoehnleProfessional GmbH+Co. KG
SOEHNLE SAS